

SYBR® Green I concentrations for real-time PCR

SL Prevost, NA Arbour, CJ Harder
Spartan Bioscience Inc.

The Spartan DX™ performs best with SYBR Green I concentrations ranging from 0.2 - 0.7 X. The optimal concentration is 0.4 X.

Introduction

SYBR Green I is a DNA-intercalating dye that is widely used for real-time PCR applications (Ref 1). It binds within the minor groove of double-stranded DNA and emits a fluorescent signal with a peak at 525 nm. Since SYBR Green I preferentially binds to double-stranded and not single-stranded DNA, it can be used to monitor the increase in double-stranded products during PCR.

There are advantages and disadvantages to using SYBR Green I for real-time PCR detection. The advantages include high detection sensitivity and low reaction cost compared to fluorescent probes. The disadvantage is that it binds to any double-stranded DNA, including non-specific primer-dimers. In addition, SYBR Green I is known to inhibit PCR at higher concentrations (Ref 2). Conversely, too low a concentration of SYBR Green I may make it difficult to detect the increase in PCR products.

The purpose of this study was to determine a useful concentration range for SYBR Green I for the Spartan DX personal DNA analyzer.

Materials and Methods

DNA extraction

DNA was extracted from clinical isolates of *Staphylococcus aureus*. For each isolate, 4-5 medium-sized bacterial colonies were re-suspended in 100 µl of lysis buffer (50 mM Tris-HCl, 50 mM NaCl, 5 mM EDTA, pH 8) with 2 µl of 1 mg/ml lysostaphin (Sigma-Aldrich, Cat No. I7386). The samples were incubated at 37°C for 30 min. Following incubation, 5 µl of 20 mg/µl Proteinase K (Sigma-Aldrich, Cat. No. p2308) was added to the mixture and tubes were shaken at 50°C for 1 h. The tubes were then incubated at 100°C for 10 min to inactivate the Proteinase K. Extracts were adjusted to 750 ng/µl, and stored at -20°C.

Real-time PCR

Oligonucleotide primers were designed against a conserved region of the bacterial *16S* gene. The forward primer sequence was 5'-cga aag cgt ggg gat caa ac-3' and the reverse primer was 5'-ccc agg cgg agt gct taa tg -3'. The primers were designed with Primer3 software (http://frodo.wi.mit.edu/cgi-bin/primer3/primer3_www.cgi). The expected amplicon

size was 125 bp.

Components of the real-time PCR amplification mixture are listed in Table 1. A 10,000 X stock of SYBR Green I (Invitrogen, Cat. No. 57563) was serially diluted so that final reaction concentrations ranged from 0.1 to 1.2 X. Samples were loaded into 0.2 ml thin-wall, flat-cap PCR tubes (Axygen, Cat. No. PCR 02C), and topped with 15 µl of mineral oil (Biotools, Cat. No. 20.032) to prevent evaporation. Real-time PCR was performed using the Spartan DX device. Table 2 shows the cycling parameters for a two-temperature program, which combines the PCR annealing and extension steps. All reactions were performed in triplicate.

DNA analysis

Fluorescence data from the real-time runs was downloaded from the Spartan DX to a computer and graphed using Microsoft Excel®. Results were confirmed by gel electrophoresis using 10 µl of the amplification products.

Results

Threshold cycles (Ct) for real-time PCR ranged from 11 to 12 with SYBR Green I concentrations of 0.2 to 0.7 X (Fig 1). Concentrations greater than 0.7 X resulted in a delay of the Ct values from 17 to 31 (Fig 2). Concentrations less than 0.2 X were difficult to distinguish from baseline fluorescence. Gel electrophoresis confirmed amplification of the expected

Component	Final amount
10X PCR Reaction Buffer (No MgCl ₂) (Invitrogen)	1 X
MgCl ₂ (Invitrogen)	2.5 mM
dNTP mix (Invitrogen)	0.125 mM
Taq DNA polymerase (Invitrogen)	1 U
SYBR Green I (Invitrogen)	0.1 - 1.2 X
PCR primers (Sigma-Aldrich)	0.5 µM
Template DNA	1.5 µg
Sterile water	
Total reaction volume	20 µl

Table 1. Components of amplification mixture.

Step	Temperature	Time	Cycles
Initial denaturation	93.6°C	30 s	1
Denaturation	93.6°C	30 s	30
Annealing/extension	56.9°C	30 s	30

Table 2. Cycling parameters.

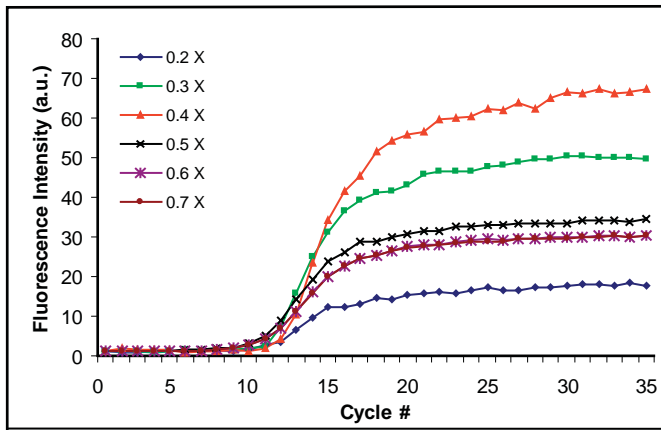


Figure 1. Real-time PCR results with SYBR Green I concentration ranging from 0.2 to 0.7 X.

125 bp fragment for samples containing 0.1 X to 0.8 X SYBR Green I (data not shown). We also observed variations in SYBR Green I fluorescence depending on the lot used.

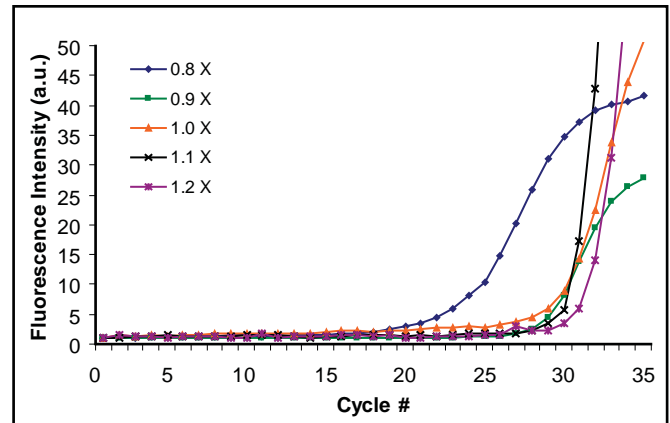


Figure 2. Real-time PCR results for SYBR Green I concentrations ranging from 0.8 to 1.2 X (delayed Ct).

Discussion and Conclusions

We observed that concentrations of SYBR Green I above 0.7 X were inhibitory to PCR as evidenced by the delay in Ct value. This was expected based on previous reports in the literature (Ref 3). Concentrations less than 0.2 X were too low to be detectable above baseline fluorescence. Therefore, a SYBR Green I concentration of 0.2 to 0.7 X appears to be a workable range for the Spartan DX. To avoid problems stemming from variations between lots, we recommend using 0.4 X SYBR Green I for most real-time PCR applications.

References

1. Monis PT, et al (2004). Comparison of SYTO9 and SYBR Green I for real-time Polymerase Chain Reaction and investigation of the effect of dye concentration on amplification and DNA melting curve analysis. *Analytical Biochemistry*. 340:24-34.
2. Epsy MJ, et al (2006). Real-Time PCR in clinical microbiology: applications for routine laboratory testing. *Clinical Microbiology Reviews*. 19(1):165-256.
3. Wittwer CT, et al (1997). Continuous fluorescence monitoring of rapid cycle DNA amplification. *Biotechniques*. 22:130-131, 134-138.

Disclaimer

PCR and real-time PCR processes are covered by patents issued and applicable in certain countries. This product is not licensed under these patents. Spartan does not encourage or support the unauthorized or unlicensed use of PCR or real-time PCR processes. Use of this instrument is recommended for persons that either have licenses to perform PCR and real-time PCR, or are not required to obtain licenses. Users interested in obtaining a license for these patents should contact the respective patent and license owners.

This product is not licensed under U.S. Patent Nos. 6,174,670 and 6,658,627, for use of SYBR Green® I in PCR. Users interested in obtaining a license for these patents should contact Idaho Technology, 390 Wakara Way, Salt Lake City, UT 84108, 801-736-6354, it@idahotech.com.

Trademarks

Spartan DX™ is a registered trademark of Spartan Bioscience Inc. All other trademarks are the sole property of their respective owners.